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
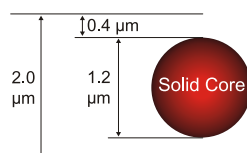
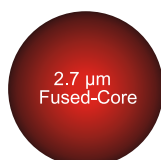
Best Fused-Core UHPLC Column
An optimized solution for high-throughput small molecule analysis

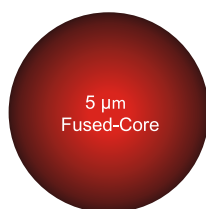Fast HPLC on ANY System
A practical solution that delivers UHPLC performance from any HPLC

The Lab Work-Horse Column
True plug and play solution for improving existing 3 or 5 µm porous particle HPLC columns

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
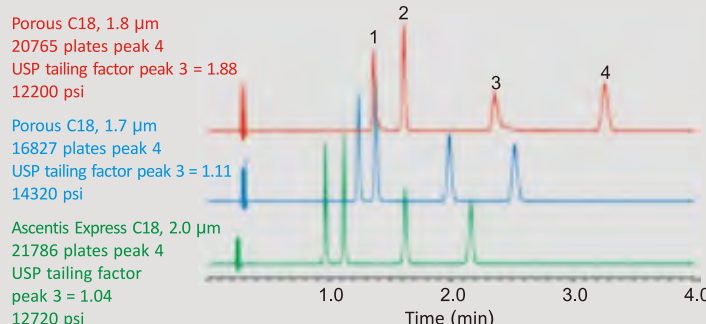
column: Ascentis Express C18, 10 cm x 2.1 mm I.D., 2.0 µm (50813-U)
mobile phase: (A) water with 0.1% formic acid; (B) acetonitrile with 0.1% formic acid; (65:35, A:B)
flow rate: 0.8 mL/min
temp.: 35 °C
det: UV, 240 nm
injection: 0.5 µL
sample: 50 µg/mL in 80:20, water:methanol
instrument: Dionex Ultimate 3000

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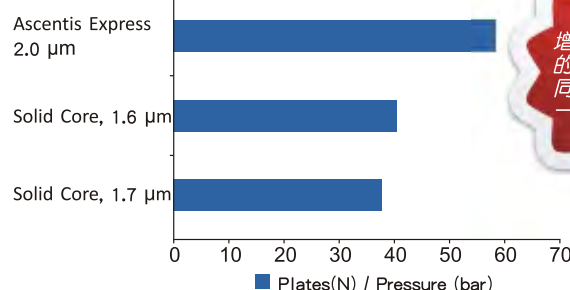Selecting Your Ascentis® Express 2.0 µm Column

Which column I.D. is best for my needs?

- If you are doing Mass Spec.....2.1 mm I.D.
- If you want high loading3.0 mm I.D.
- If you want high sensitivity.....2.1mm I.D.
- If you want high performance.....3.0 mm I.D.

Which column length is best for my needs?

- If you want to maximize the speed of your application2 to 5 cm
- If you want a balance of resolution and speed10 cm

What flow rate is best for my needs?

- If you are using a column with a 3.0 mm I.D.....0.8 to 2.0 mL/min*
- If you are using a column with a 2.1 mm I.D.....0.4 to 1.8 mL/min*

*Higher flow rates may be allowed by certain instruments.

Ascentis Express 2.0 µm Fused-Core Silica Characteristics 粒子規格

Pore Size	90Å
Max Temp.....	60 °C
Pressure	1000bar (14500 psi)
Features	Endcapped
Operating pH Range.....	2-9
Surface Area.....	120 m ² /g

Get Started

Additional resources are available for helping you implement Ascentis Express HPLC Columns into your laboratory.

For product information,webinars, ordering and real time availability information, visit
sigma-aldrich.com/express

Our technical service staff is ready to answer questions.
EU: eurtechserv@sial.com
US: techserv@sial.com

A technical seminar can be arranged on-site or via the web.
Request viaseminars@sial.com

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Description	Pkg. Size	Cat. No.
Universal Guard Holder		
Holder w/EXP Titanium Hybrid Ferrule (cartridge not included)	1	53500-U

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UNI-ONWARD Corp.