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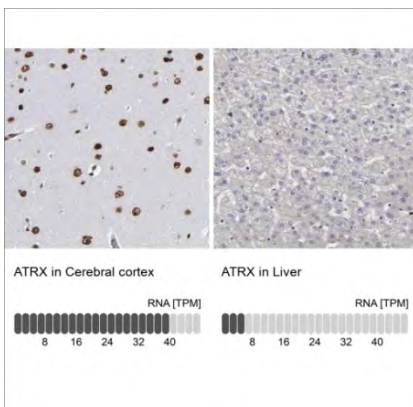
Fig1: orthogonal validation of [Anti-ATRX antibody \(HPA001906\)](#) in IHC by staining of cerebral cortex and liver tissues, comparing to ATRX RNA-seq data are presented for the same tissues.

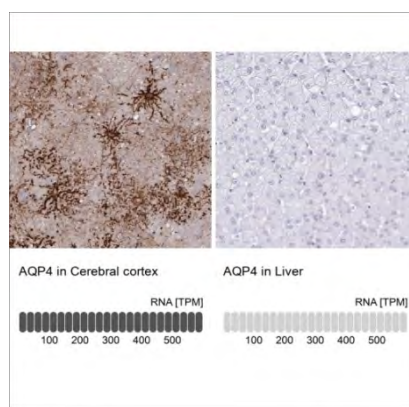
Fig2: orthogonal validation of [Anti-AQP4 antibody \(HPA014784\)](#) in IHC by staining of cerebral cortex and liver tissues, comparing to ATRX RNA-seq data are presented for the same tissues.

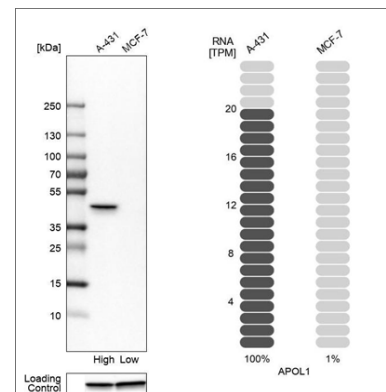
Fig3: orthogonal validation of [Anti-APOL1 antibody \(HPA018885\)](#) by Western blot analysis in human cell lines A-431 and MCF-7. Corresponding APOL1 RNA-seq data are presented for the same cell lines

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

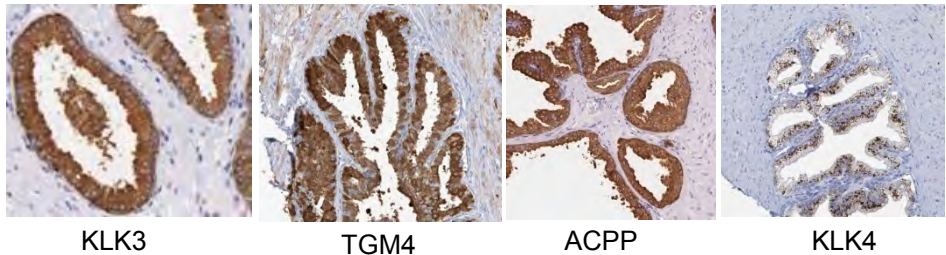
ACPP

KLK4

Target protein	品號	應用	Description	Predicted localization
KLK3	HPA000764	IHC	kallikrein-related peptidase 3	Secreted
TGM4	HPA032072	IHC	transglutaminase 4	Intracellular
ACPP	HPA063916	IHC, WB	acid phosphatase, prostate	Intracellular,Membrane, Secreted
KLK4	HPA051839	IHC	kallikrein-related peptidase 4	Intracellular,Secreted
SLC45A3	HPA019075	IHC,IF,WB	solute carrier family 45, member 3	Membrane
MSMB	HPA051257	IHC, WB	microseminoprotein, beta-	Secreted

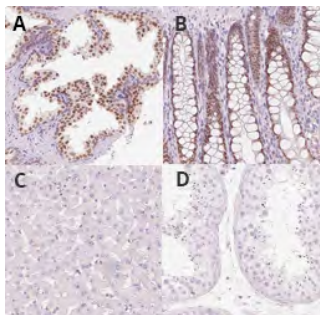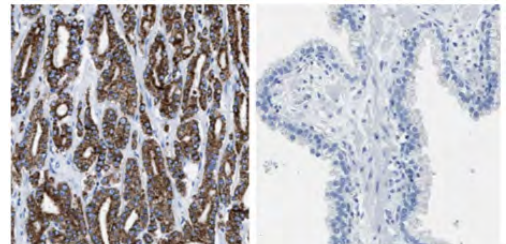IHC staining of human prostate (A), colon (B), liver (C) and testis (D) tissues with [Anti-HOXB13 antibody \(HPA062852\)](#) showing grouped expression in prostate and colon.

The [Anti-AMACR antibody \(HPA019527\)](#) shows strong positivity in prostate cancer samples (left) compared to normal prostate (right).

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
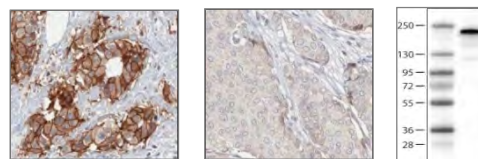
HER2 /ERBB2


Immunohistochemical staining of human breast tumour using Anti-HER2 (AMAb90627) shows strong membranous (combined with moderate cytoplasmic) positivity in tumour cells in HER2-positive ductal carcinoma, while HER2-negative ductal carcinoma shows no membranous positivity. By WB analysis, HER2 is detected in the breast cancer cell line SK-BR-3.

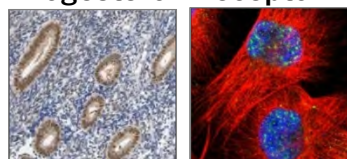Progesteron receptor


IHC staining using the Anti-PGR antibody (HPA004751) in normal human corpus (uterine) tissue shows strong nuclear positivity in glandular cells. In the presented breast cancer sample, the staining of tumor cells is also nuclear. ICC-IF shows nuclear staining in U-251MG cells.

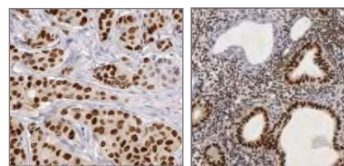Estrogen Receptor


IHC staining using the Anti-ESR1 antibody (HPA000450) shows strong nuclear positivity in glandular and stromal cells of human corpus, uterine tissue and in tumor cells in breast cancer.

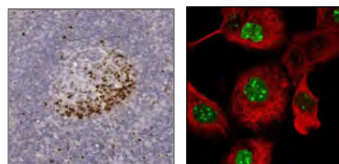Ki67


IHC staining of human tonsil tissue using the Anti-MKI67 antibody (HPA001164) shows nuclear staining of reaction center cells

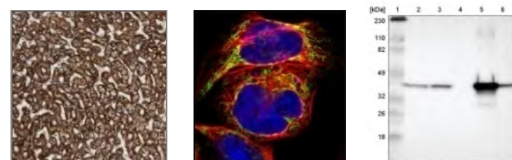ACAT1 (HPA004428)

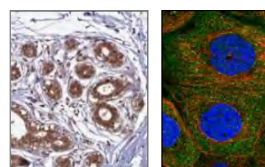
Immunoperoxidase staining of formalin-fixed, paraffin-embedded human duodenum tissue showing granular cytoplasmic staining of glandular cells.

BRCA1

The Anti-BRCA1 antibody (HPA034966) shows positivity in glandular cells in normal human breast tissue and in tumor cells in breast cancer samples using IHC.

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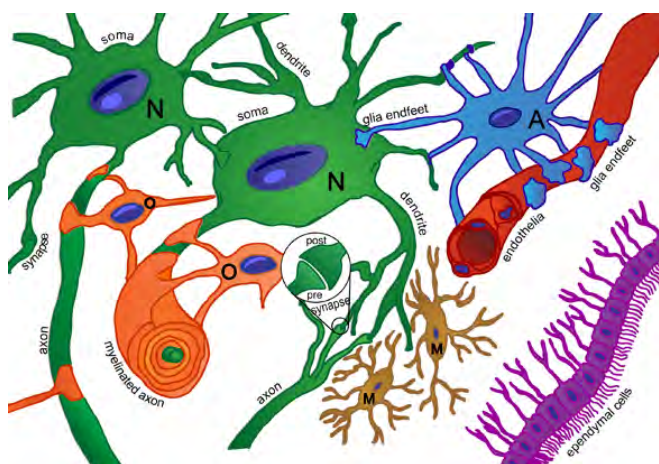The different types of neurological cell classes (ependymal cells, purple), subpopulations (A=astrocytes, blue; M=microglia, brown; N=neuron, green; O=oligodendrocytes, orange), and subcellular locations (axon, dendrite, synapse, and glia endfeet).

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